


Analysis of 13 kinds of Aldehydes (Detected by Pre-Column Method with DNPH)

Data No. LA960-0894

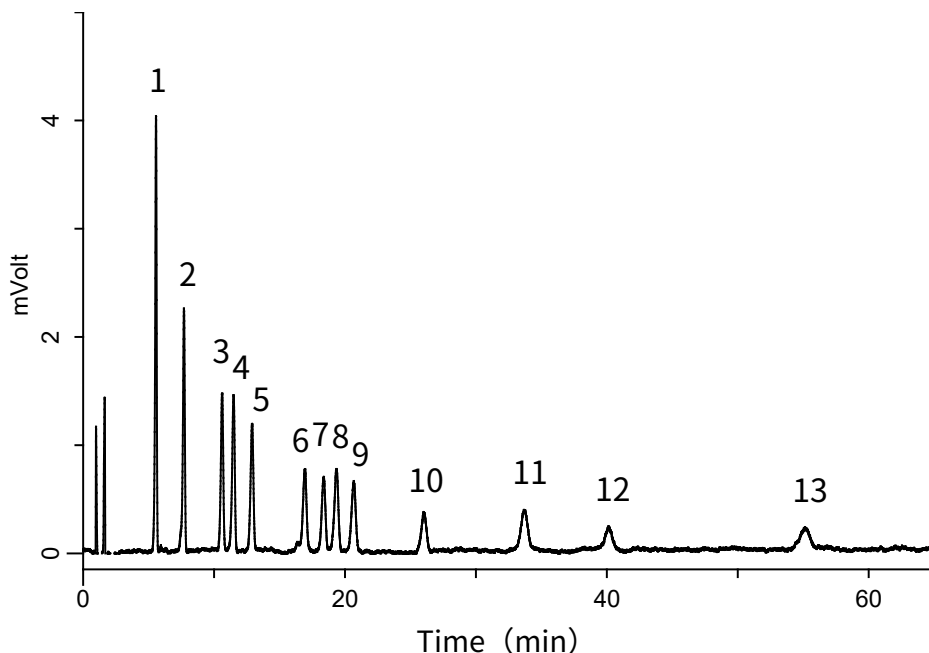

Conditions

System : LC800 HPLC system
Column : InertSustain C18 HP (3μm, 250 x 3.0 mm I.D.)
Column Cat. No. : 5020-14426
Eluent : A) CH₃CN
 : B) H₂O
 : C) THF
 : A/B/C = 35/55/10, v/v/v
Flow rate : 1.0 mL/min
Col. Temp. : 40 °C
Detection : UV 360 nm (LC800 UV Detector)
Injection Vol. : 10 μL

Analyte:

1. DNPH-Formaldehyde	150 μg/L
2. DNPH-Acetaldehyde	150 μg/L
3. DNPH-Acetone	150 μg/L
4. DNPH-Acrolein	150 μg/L
5. DNPH-Propionaldehyde	150 μg/L
6. DNPH-Crotonaldehyde	150 μg/L
7. DNPH-Methylethylketone	150 μg/L
8. DNPH-Methacrolein	150 μg/L
9. DNPH- <i>n</i> -Butyraldehyde	150 μg/L
10. DNPH-Benzaldehyde	150 μg/L
11. DNPH- <i>n</i> -Valeraldehyde	150 μg/L
12. DNPH- <i>m</i> -Tolualdehyde	150 μg/L
13. DNPH-Hexanal	150 μg/L