
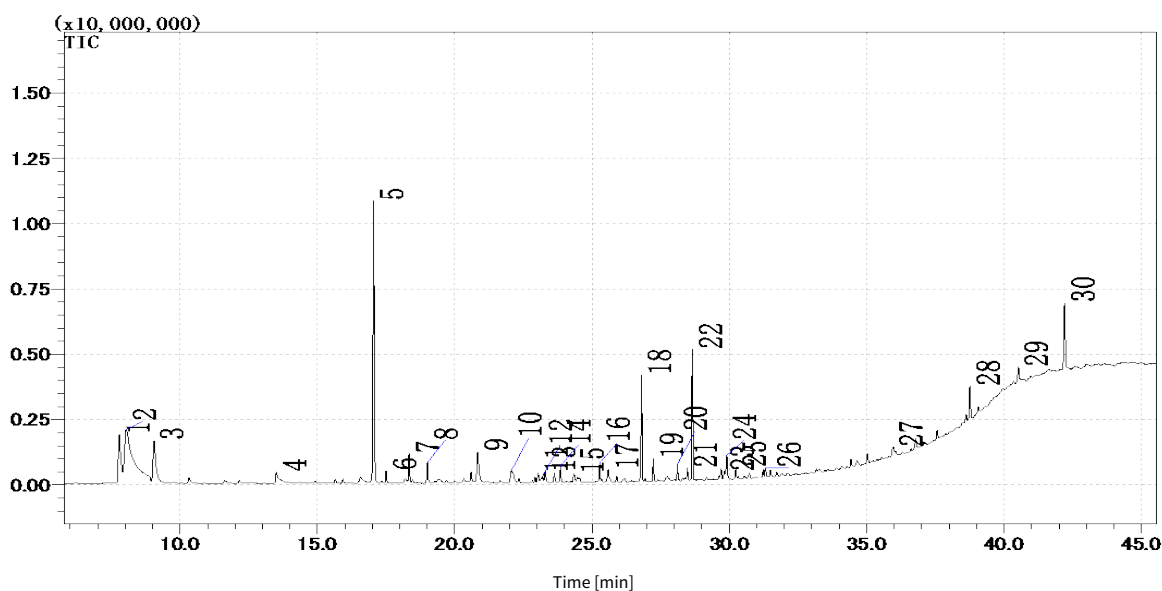


Handy TD and MonoTrap RGC18TD were used to screen the flavor components of Mikan juice. Handy TD is a compact, portable and easy-to-operate thermal desorption device. After sample collection, place the MonoTrap in the glass liner in the Handy TD, and introduce the heat-desorbed components into the GC inlet. A large number of terpenes, including Limonene, were found in citrus fruit.

Preliminary processing procedure

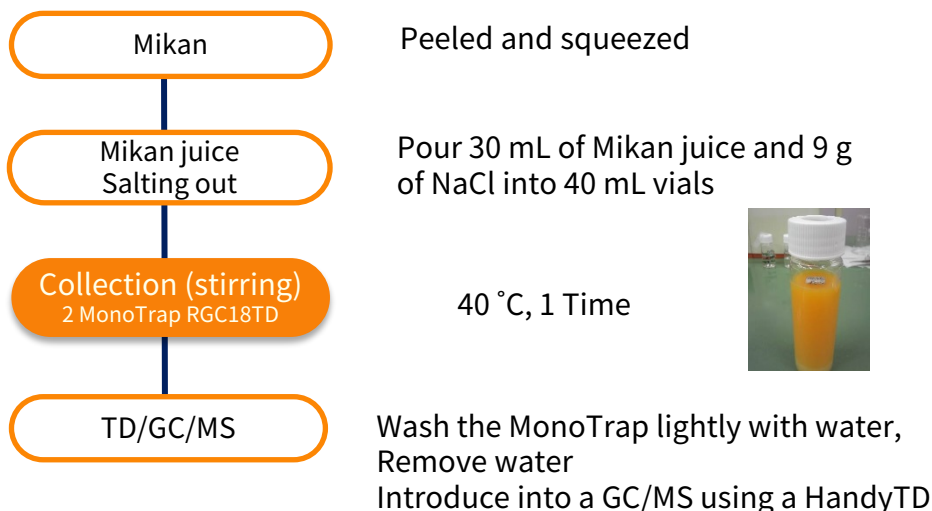

* Standard samples are not used for qualitative analysis.
Results from a library search.

- | | | |
|-------------------|--------------------|---------------------------|
| 1. Ethyl Acetate | 11. Acetic acid | 21. Citral |
| 2. Methyl Alcohol | 12. Linalool oxide | 22. Terpeneol |
| 3. Ethanol | 13. Heptanol | 23. Isopiperitenol |
| 4. Hexanal | 14. Linalool oxide | 24. Isopiperitenol |
| 5. D-Limonene | 15. Formic acid | 25. Dibutyl formamide |
| 6. 2-Hexenal | 16. Linalool | 26. Carveol |
| 7. Terpinene | 17. Octanol | 27. Dihydroxyacetone |
| 8. Cymene | 18. Terpeneol | 28. Dihydroxy maltol |
| 9. 1-Hexanol | 19. Menthadienol | 29. Dihydroactinidiolide |
| 10. Nonanal | 20. Menthadienol | 30. Hydroxymethylfurfural |

Product used

Portable thermal desorber HandyTD TD265

Product name : HandyTD TD 265
Cat.No. : 2709-80000



Collection MonoTrap RGC18TD



MonoTrap RGC18TD

Product name : MonoTrap RGC 18 TD
Cat.No. : 1050-74201



MonoTrap Trial Kit for TD

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome
Shinjuku-ku, Tokyo
163-1130, Japan

Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
USA

Phone: +1-310-265-4424
Fax: +1-310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences B.V.

Dillenburgstraat 7C
5652AM, Eindhoven
The Netherlands

Phone: +31-40-254-9531
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (Shanghai) Limited

Tower B, Room 2003
Far East International Plaza
No.317 Xianxia Road, Changning District
Shanghai, China 200051

Phone: +86-21-62782272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn



GL Sciences
International Distributors

Visit our Website at www.glsciences.com/distributors