

Concentration Analysis of Volatile Components in Wasabi Paste - Using HandyTD Portable Thermal Desorber and MonoTrap Simplified Enrichment Tools

The volatile components of commercial GPS TD. HandyTD is a compact, portable and easy-to-operate thermal desorption device. After sample wasabi paste were screened and analyzed using HandyTD and a simplified enrichment tool MonoTrap Rollection, set a glass liner containing the MonoTrap and introduce the heat-desorbed components into the GC inlet. A large number of different characteristic sulfur compounds were detected.

Pretreatment procedure

Wasabi paste

Place 6.25 g in a 44-mL vial

Collection (HS)
One MonoTrap RGPS TD

37 °C for 1 hour




TD/GC/MS

GC Conditions

System : GC - MS - Thermal Desorption (HandyTD TD265)
Column : InertCap Pure-WAX 0.32 mm I.D. x 60 m, df = 0.50 µm
Col.Cat. No. : 1010-68264
Col.Temp. : 40 °C (5 min) - 4 °C/min - 250 °C
Carrier Gas : He, 1.88 mL/min
GC Inlet : 250 °C, Split 7:1
Detection : MS Scan (*m/z* 15-450)

HandyTD Conditions

Desorb Temp. : 40 °C - 45 °C/sec - 250 °C (1.5 min)
Pre Desorb Press. : 108 kPa

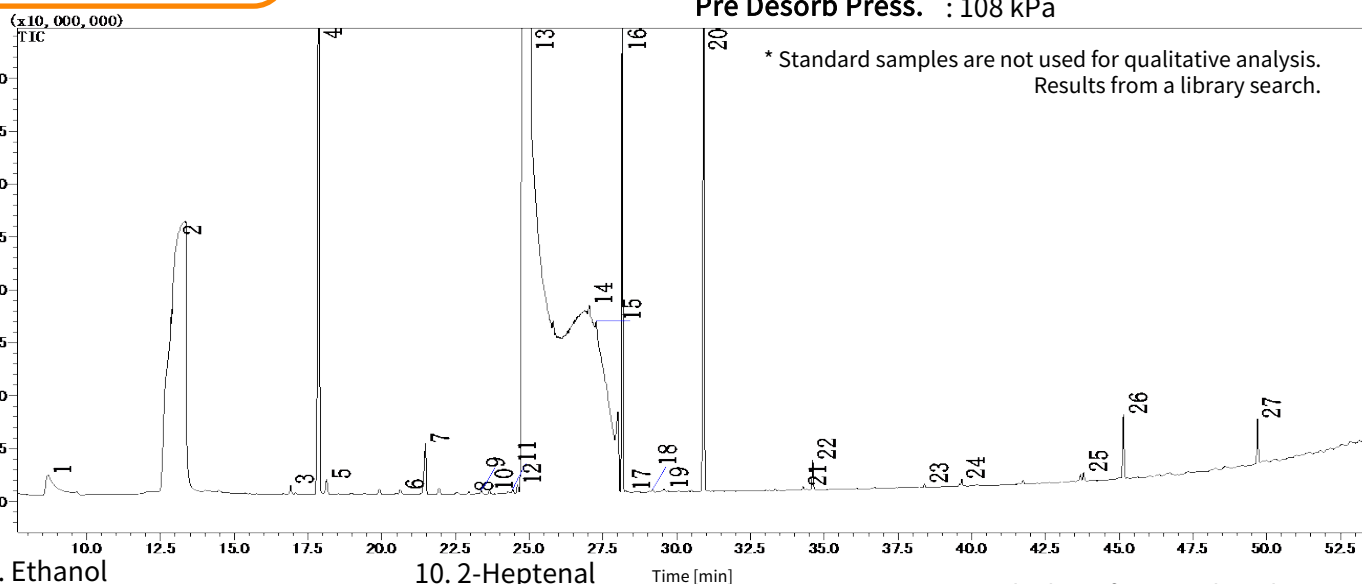

- | | | |
|--------------------------------|--------------------------------------|----------------------------------|
| 1. Ethanol | 10. 2-Heptenal | 19. Tetrahydro-2-furanmethanol |
| 2. Water | 11. 5-Cyano-1-pentene | 20. Unknown |
| 3. 2-Butenenitrile | 12. 1-Hexanol | 21. Dihydro-2(3H)-thiophenone |
| 4. 1-Pentyne | 13. Allyl isothiocyanate | 22. Unknown |
| 5. Isopropyl isothiocyanate | 14. 2-Methylbutyl isothiocyanate | 23. 4-(Methylthio) butanenitrile |
| 6. 1-Pentanol | 15. 1-Isothiocyanato-3-methyl-butane | 24. Hexanoic acid |
| 7. 2-Isothiocyanatbutane | 16. 4-Isothiocyanato-1-butene | 25. Iberverin |
| 8. 1-(1-Ethoxyethoxy)-3-hexene | 17. Furfural | 26. Benzenepropanenitrile |
| 9. Isobutyl isothiocyanate | 18. Pentyl isothiocyanate | 27. Phenylethyl isothiocyanate |

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome
Shinjuku-ku, Tokyo
163-1130, Japan

Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
USA

Phone: +1-310-265-4424
Fax: +1-310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences B.V.

Dillenburgstraat 7C
5652AM, Eindhoven
The Netherlands

Phone: +31-40-254-9531
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (Shanghai) Limited

Tower B, Room 2003
Far East International Plaza
No.317 Xianxia Road, Changning District
Shanghai, China 200051

Phone: +86-21-62782272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn



International Distributors

Visit our Website at www.glsciences.com/distributors