


Analysis of Terbinafine Chloride cream (Under the Condition of Japanese Pharmacopoeia)

Data No. LA941-811

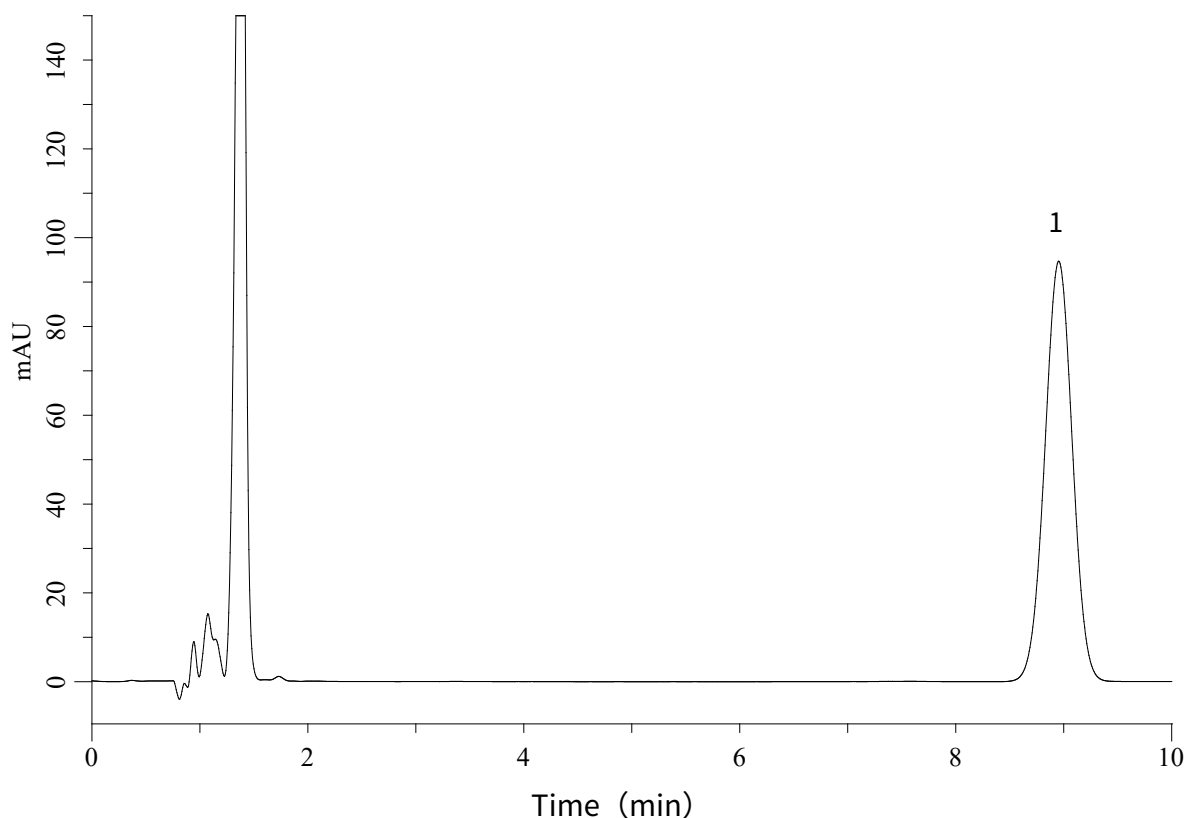

Conditions

System : GL-7400 HPLC system
Column : Inertsil WP300 C18 (5 μ m, 125 x 4.0 mm I.D.)
Column Cat. No. : 5020-87037
Eluent : A) CH₃CN
 B) THF
 C) Solition*
 A/B = 2/1/2, v/v/v

Flow rate : 1.3 mL/min
Col. Temp. : 25 °C
Detection : UV 282 nm (GL-7452A PDA Detector)
Injection Vol. : 10 μ L

Sample : Terbinafine ointment

*Dissolve 9 mL of Tetramethylammonium hydroxide (10 % in H₂O) in 2000 mL of water.
Adjust to pH 8.0 with phosphoric acid (1 in 25).

Analyte:

1. Terbinafine Chloride