

Simple Concentration and Analysis of Volatile Compounds in Socks after Wearing – Using Sorptive Media MonoTrap

A simple concentration assay of volatile compounds in socks immediately after wearing the product was carried out using MonoTrap RGC 18TD a simple concentration tool. The MonoTrap RGC18TD is packaged individually in ampoules, no pre-conditioning is required and there is no risk of contamination.

Preliminary processing procedure

Sock

Men in their 40 s worn daily

Collection (passive)
2 MonoTrap RGC18TD


In Tedler bags 37 °C for 1 hour



TD-GC-MS

Conditions

System	: GC-MS-Thermal Desorption
Column	: InertCap Pure-WAX 0.25 mm I.D x 30 m df = 0.25 µm
Col.Temp.	: 40 °C (5 min) - 10 °C/min - 250 °C
Carrier Gas	: He 1 mL/min (constant flow)
Desorb Temp.	: 200 °C
Time	: 5 min
Flow	: 2 mL/min
Split	: Splitless
Cryo Trapping	: -120 °C
Injection Temp.	: 250 °C
Detection	: MS Scan (<i>m/z</i> : 30 - 600)

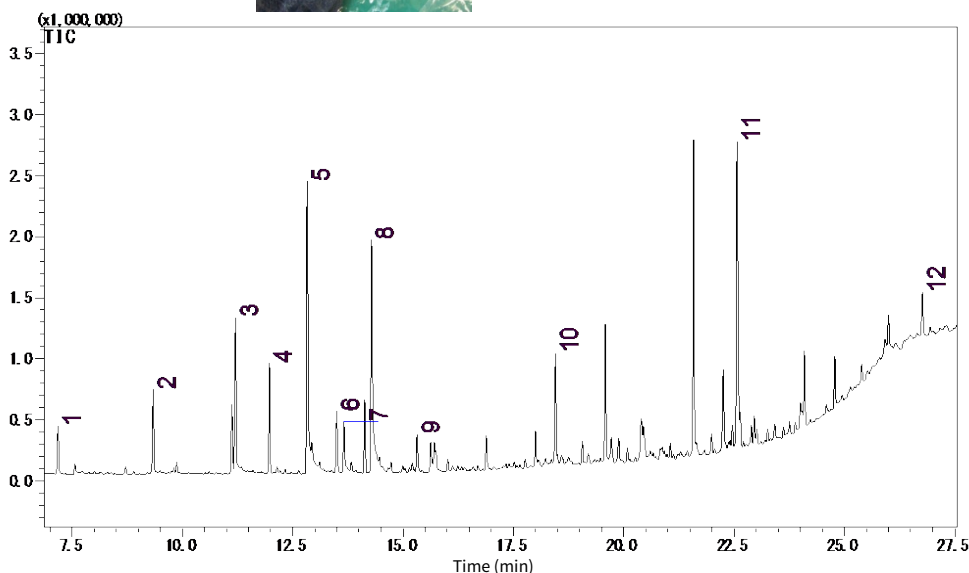

- | | |
|-----------------------------|----------------------------------|
| 1. Hexanal | 7. Dodecanal |
| 2. Heptanal | 8. Decanal |
| 3. Octanal | 9. Propylene Glycol |
| 4. 6-Methyl-5-heptene-2-one | 10. <i>Trans</i> -Geranylacetone |
| 5. Nonanal | 11. Muskalactone |
| 6. Dichlorobenzene | 12. Benzyl salicilate |

* Standard samples are not used for qualitative analysis.
Results from a library search.

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome
Shinjuku-ku, Tokyo
163-1130, Japan

Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsciences.co.jp
Web: www.glsciences.com

GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
USA

Phone: +1-310-265-4424
Fax: +1-310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences B.V.

Dillenburgstraat 7C
5652AM, Eindhoven
The Netherlands

Phone: +31-40-254-9531
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (Shanghai) Limited

Tower B, Room 2003
Far East International Plaza
No.317 Xianxia Road, Changning District
Shanghai, China 200051

Phone: +86-21-62782272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn



International Distributors

Visit our Website at www.glsciences.com/distributors