
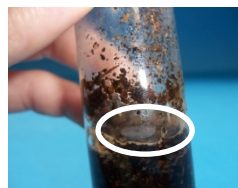


Simple Concentration Analysis of Pu-erh Tea Flavor Components - Using Sorptive Media MonoTrap

MonoTrap DCC18 (with activated carbon) was used to simply concentrate and analyze the fragrance components in pu-erh tea. Caffeine, a compound with high-boiling point was detected with high sensitivity by collecting tea leaves extracted with hot water directly on a MonoTrap sampler.

Preliminary processing procedure

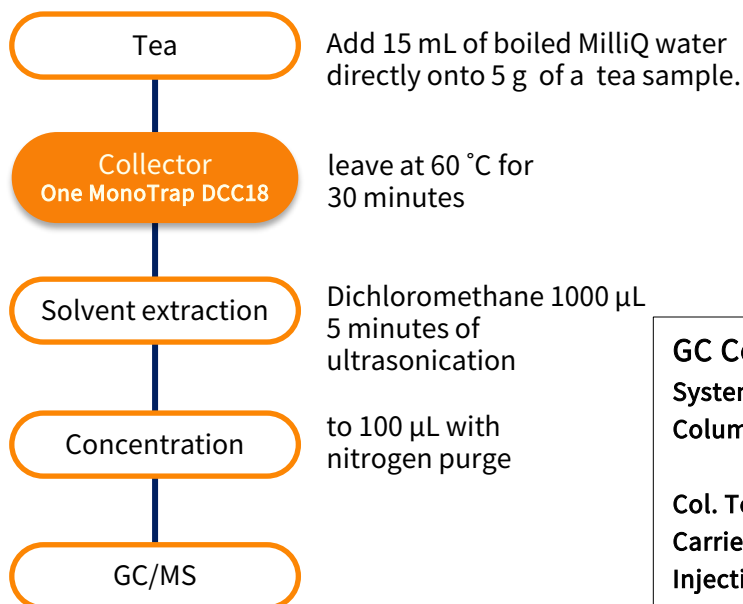


GC Conditions

System : GC - MS
Column : InertCap Pure-WAX 0.25 mm I.D. x 30 m
 df = 0.25 µm
Col. Temp. : 40 °C (5 min) - 4 °C /min - 250 °C
Carrier Gas : He 1 mL/min
Injection : Splitless
 250 °C
Detection : MS Scan (*m/z* 40-600)
Sample Size : 1 µL

MonoTrap Series

Cat.No	Description	Configuration	Size	Activated Carbon	Functional Group
1050-72101	MonoTrap DCC18	Disk	O.D.10mmx thick 1 mm	Yes	C18
1050-72201	MonoTrap RCC18	Rod	O.D.2.9 mmx I.D.1 mm x Height 5 mm	Yes	C18
1050-71101	MonoTrap DSC18	Disk	O.D.10 mm x thick 1 mm	No	C18
1050-71201	MonoTrap RSC18	Rod	O.D.2.9 mm x I.D.1 mm x Height 5 mm	No	C18

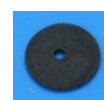

DCC18



RCC18




DSC18




RSC18

Initial use of the convenient MonoTrap and overview of the start-up kit

	Description	Qty.
①	MT Holder	5
②	MT Stand	1
③	MT Extract Cup with Vial(20 mL)	5
④	Clean Pin Hole Septum with vial(40 mL)	5
⑤	200 µL glass insert (flat bottom)	40
⑥	MonoTrap DCC18	20
⑦	MonoTrap RCC18	20
⑧	MonoTrap DSC18	20
⑨	MonoTrap RSC18	20

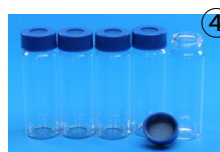

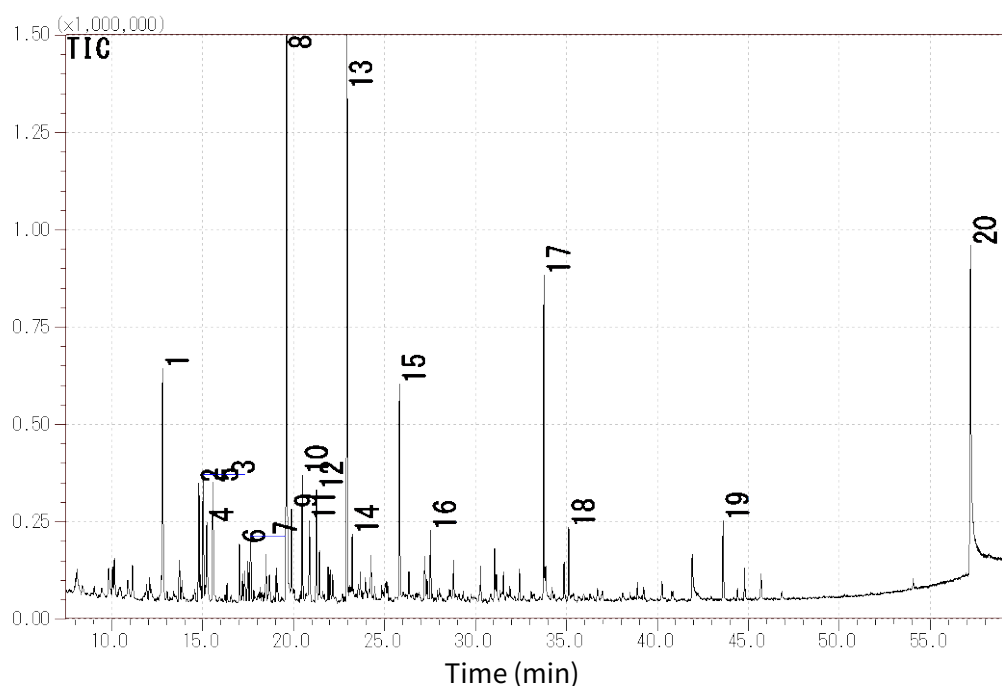

- | | |
|-----------------------------|----------------------------|
| 1. Methylpyrazine | 11. 2-Acetylfuran |
| 2,3. Dimethylpyrazine | 12. Benzaldehyde |
| 4. Ethylpyrazine | 13. 2-Formyl-5-methylfuran |
| 5. 6-Methyl-5-heptene-2-one | 14. Methyl 2-furoate |
| 6. 2-Ethyl-6-methylpyrazine | 15. 2-Furanmethanol |
| 7. Trimethylpyrazine | 16. Dimethoxybenzene |
| 8. Furfural | 17. Trimethoxybenzene |
| 9. Acetol acetate | 18. Pyrrole-2-aldehyde |
| 10. 2,4-Heptadien-1-al | 19. Coumaran |
| | 20. Caffein |

* According to library search results

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome
Shinjuku-ku, Tokyo
163-1130, Japan

Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
USA

Phone: +1-310-265-4424
Fax: +1-310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences B.V.

Dillenburgstraat 7C
5652AM, Eindhoven
The Netherlands

Phone: +31-40-254-9531
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (Shanghai) Limited

Tower B, Room 2003
Far East International Plaza
No.317 Xianxia Road, Changning District
Shanghai, China 200051

Phone: +86-21-62782272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn



International Distributors

Visit our Website at www.glsciences.com/distributors