

Concentration analysis of volatile constituents of mosquito incense coil - Using Handy TD Portable Thermal Desorber and MonoTrap Simple Enrichment Tools

Handy TD and MonoTrap RGPSTD were used to screen the volatile components from the aroma of a mosquito-coil. Handy TD is a compact, portable, and easy-to-operate thermal desorption device. After collection, place the MonoTrap in a glass liner in the Handy TD, and introduce the heat-desorbed components into the GC inlet. A number of pyridines and aromatic compounds were detected.

Preliminary processing procedure

Mosquito coil

3.62 g

Collection (HS)
3 MonoTrap RGPSTDAt room
temperature
for 1 hour

TD/GC/MS


Introduced into GCMS using HandyTD

GC Conditions

System : GC - MS - Handy TD TD265
Column : InertCap Pure-WAX
 0.25 mm I.D. x 60 m, df = 0.25 μ m
Col.Temp. : 40 °C (5 min) - 6 °C/min - 250 °C (10min)
Carrier Gas : He, 1.38 mL/min (149.6 kPa at 40 °C)
GC Inlet : 250 °C, Split 9:1
Detection : MS Scan (m/z : 45-450)

HandyTD Conditions

DesorbTemp. : 40 °C - 45 °C/sec-250 °C (1.5 min)
Pressure : 180 kPa

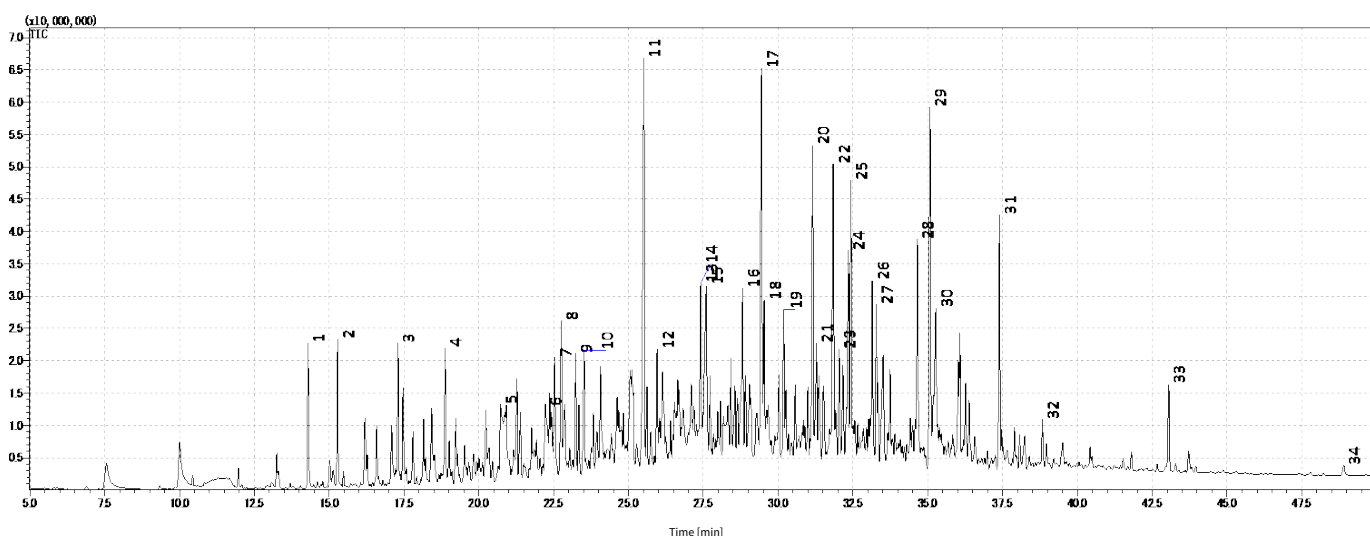

* Standard samples are not used for qualitative analysis.
Results from a library search.

- | | | |
|------------------------------|----------------------------|--------------------------|
| 1. Pyridine | 12. Methyl imidazole | 23. Dimethyl naphthalene |
| 2. Methyl pyridine | 13. Azulene | 24. Indanone |
| 3. Methyl pyridine | 14. Azulene | 25. Ethyl guaiacol |
| 4. Octanenitrile | 15. Dihydropyran | 26. Cresol |
| 5. Acetic acid | 16. Methyl cyclopentadione | 27. Cresol |
| 6. 2,3-Butanediol, dinitrate | 17. Ethyl furaldehyde | 28. Eugenol |
| 7. Trimethyl decane | 18. Methyl naphthalene | 29. Vinyl guaiacol |
| 8. Methyl cyclopentenone | 19. Methyl naphthalene | 30. Methylimidazole |
| 9. Methyl cyclopentenone | 20. Creosol | 31. Ioeugenol |
| 10. Pentadecene | 21. Maltol | 32. Indole |
| 11. Furanmethanol | 22. Phenyl carbamate | 33. Carboxamide |
| | | 34. Hydroxy phenol |

Product used

Portable thermal desorber HandyTD TD265

Product name : HandyTD TD 265
Cat.No. : 2709-80000



Collection MonoTrap RGC18TD



MonoTrap RGC18TD

Product name : MonoTrap RGC 18 TD
Cat.No. : 1050-74201



MonoTrap Trial Kit for TD

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome
Shinjuku-ku, Tokyo
163-1130, Japan

Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
USA

Phone: +1-310-265-4424
Fax: +1-310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences B.V.

Dillenburgstraat 7C
5652AM, Eindhoven
The Netherlands

Phone: +31-40-254-9531
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (Shanghai) Limited

Tower B, Room 2003
Far East International Plaza
No.317 Xianxia Road, Changning District
Shanghai, China 200051

Phone: +86-21-62782272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn



International Distributors

Visit our Website at www.glsciences.com/distributors