

Easy Concentration & Analysis of Roasted Seaweed by MonoTrap - Thermal Desorption

Introduction and analysis of roasted seaweed by Thermal Desorption using MonoTrap for TD. Compounds such as di-methyl sulfide and β -Ionone recognized as the smell of ocean were detected.

Protocol

Roasted seaweed

Roast and cut the seaweed

Head Space Sampling
MonoTrap RGC18 TD×1pcs



At 40 °C for 3 hours

TD/GC/MS



System	: GC/MS-Thermal Desorption (T-Dex II)
Column	: InertCap Pure-WAX 0.25 mm I.D. × 30 m, df = 0.25 μ m
Col.Temp.	: 40 °C (5 min) – 6 °C/min – 250 °C
Carrier Gas	: He, 1 mL/min (constant flow)
Desorb Temp.	: 200 °C
Time	: 5 min
Flow	: 5 mL/min
Split	: Splitless
Cryo Trapping	: -150 °C
Injection Temp.	: 250 °C
Detection	: MS Scan (m/z 28.5 - 600) : SIM (62,61,47 for Dimethyl sulfide)


<Scan>

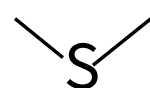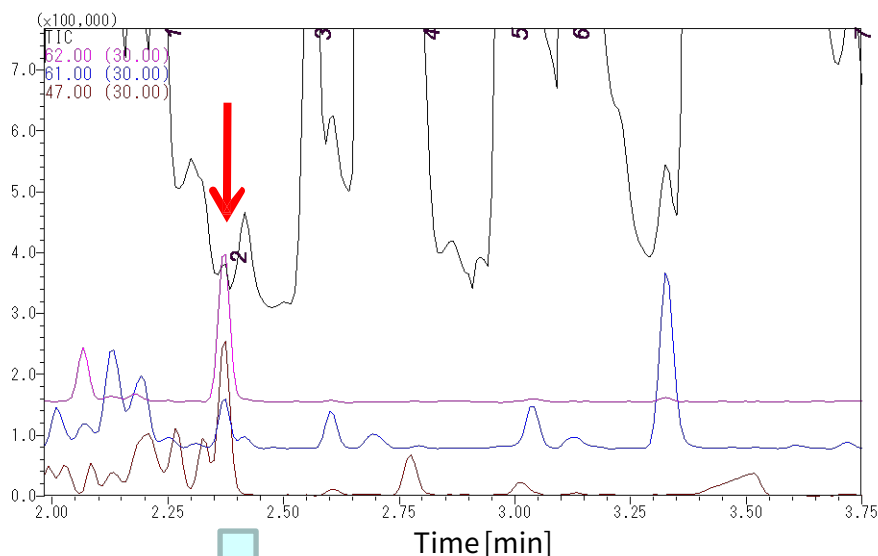


Dimethyl sulfide

m/z 62,61,47

<SIM>

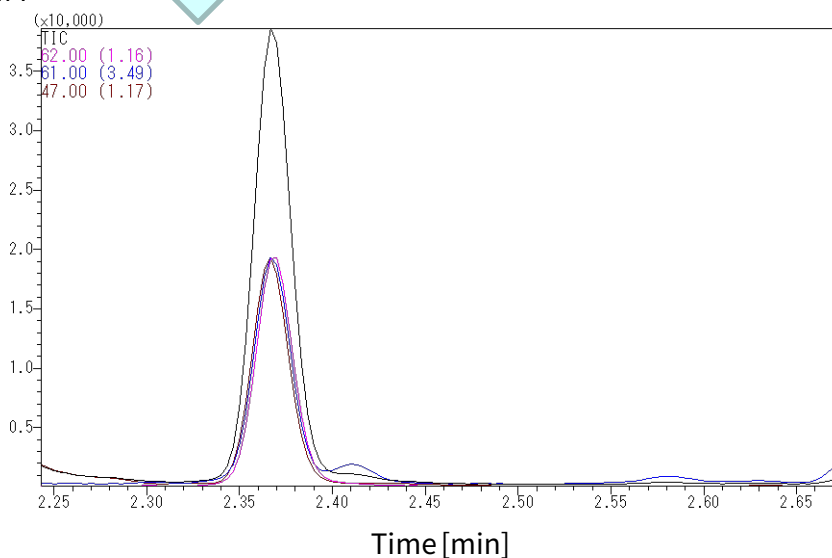

Dimethyl sulfide was
well detected by SIM

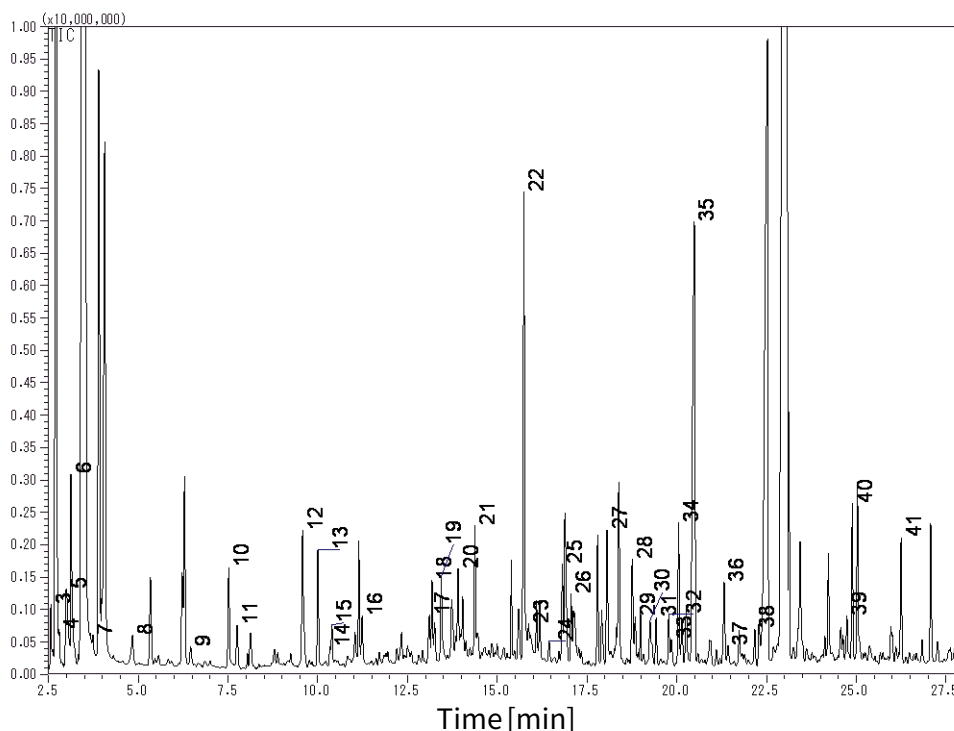

- | | |
|------------------------------------|---|
| 1. Acetaldehyde | 22. Nonanal |
| 2. Dimethyl sulfide | 23. Trimethylpyrazine |
| 3. Propanal | 24. 2-Octenal |
| 4. Methyl acetate | 25. 2-Ethyl-3,6-dimethylpyrazine |
| 5. Trimethylamine | 26. 1-Octen-3-ol |
| 6. Trimethylamine | 27. Decanal |
| 7. Isovaleraldehyde | 28. 2-Nonenal |
| 8. Pentanal | 29. 3-Caranol |
| 9. 1-Propanol | 30. Dimethyl Sulfoxide |
| 10. Hexanal | 31. 3,5-Octadien-2-one |
| 11. 2-Methyl-2-butenal | 32. 2,6-Nonadienal |
| 12. 2-Ethyl-trans-2-butenal | 33. Propylene Glycol |
| 13. 1-Penten-3-ol | 34. 2,6-Dimethylcyclohexanol |
| 14. 2-Heptanone | 35. Butyrolactone |
| 15. Heptanal | 36. N-Methyl-2-pyrrolidinone |
| 16. 2-Hexenal | 37. 2(5H)-Furanone, 5-methyl- |
| 17. Acetoin | 38. 2-Dodecanone |
| 18. Octanal | 39. α -Ionone |
| 19. Acetol | 40. <i>trans</i> -Geranylacetone |
| 20. 2,5-Dimethylpyrazine | 41. β-Ionone |
| 21. 6-Methyl-5-heptene-2-one | |

*Red: Compounds of Roasted Seaweed

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome
Shinjuku-ku, Tokyo
163-1130, Japan

Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
USA

Phone: +1-310-265-4424
Fax: +1-310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences B.V.

Dillenburgstraat 7C
5652AM, Eindhoven
The Netherlands

Phone: +31-40-254-9531
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (Shanghai) Limited

Tower B, Room 2003
Far East International Plaza
No.317 Xianxia Road, Changning District
Shanghai, China 200051

Phone: +86-21-62782272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn



International Distributors

Visit our Website at www.glsciences.com/distributors