


Analysis of Pazufloxacin mesilate

(Under the Condition of the draft for the Japanese Pharmacopoeia)

Data No. LB444-0894

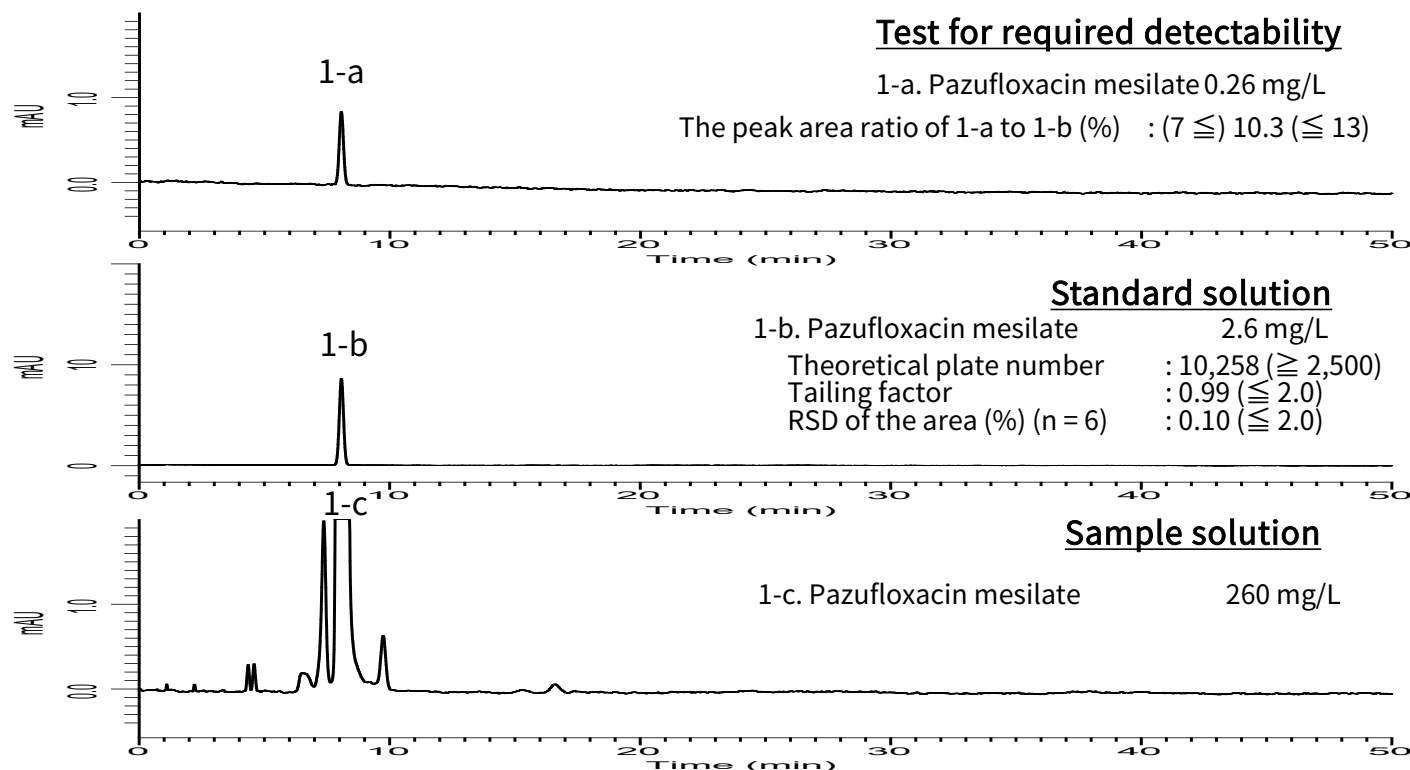

Conditions

System : GL7700 HPLC system
Column : Inertsil ODS-2
 (5 μ m, 150 x 4.6 mm I.D.)
Column Cat. No. : 5020-01124
Eluent : Buffer*
Flow Rate : 1.57 mL/min
Col. Temp. : 40 °C
Detection : UV 254 nm (PD7752 PDA Detector)
Injection Vol. : 20 μ L
Sample : Standard

Analyte:

1. Pazufloxacin mesilate

*Dissolve 1.08 g of sodium 1-octanesulfonate in 1000 mL of a mixture of diluted phosphoric acid (1 in 1000) and acetonitrile (39:11).