


Analysis of Glycyrrhiza

(Under the Condition of the Japanese Pharmacopoeia 17th edition)

Data No. LB482-0919

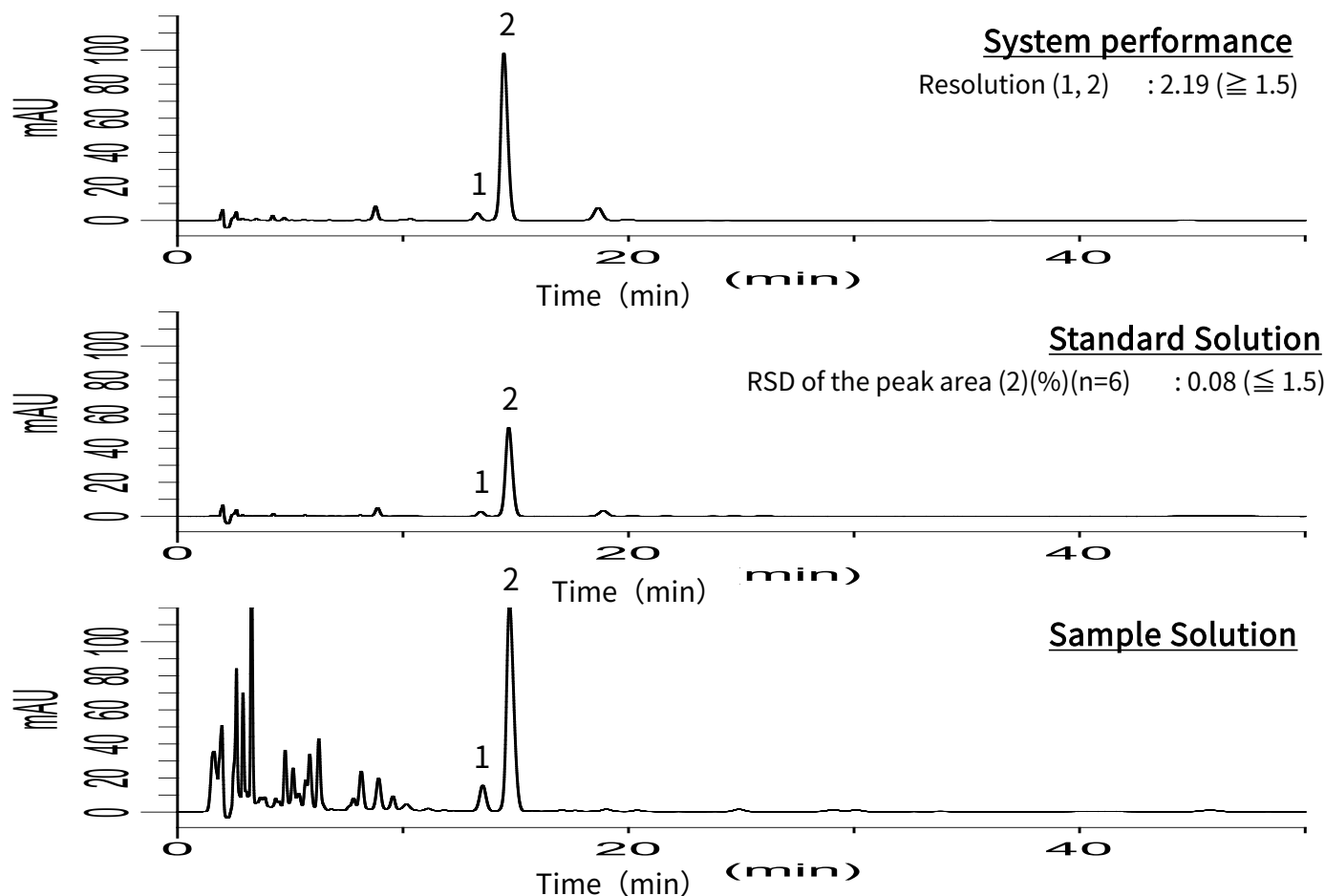

Conditions

System : GL7700 HPLC system
Column : Inertsil ODS-4
(5 μ m, 150 x 4.6 mm I.D.)

Column Cat. No. : 5020-03945

Eluent : Solution*

Flow Rate : 0.9 mL/min

Col. Temp. : 40 °C

Detection : UV 254 nm (PD7752 PDA Detector)

Injection Vol. : 10 μ L

*Dissolve 3.85 g of ammonium acetate in 720 mL of water, and add 5 mL of acetic acid and 280 mL of acetonitrile.

Analyte:

1. Impurity
2. Glycyrrhizic acid