


Analysis of Eplerenone

(Under the Condition of the draft for the Japanese Pharmacopoeia)

Data No. LB413-0894

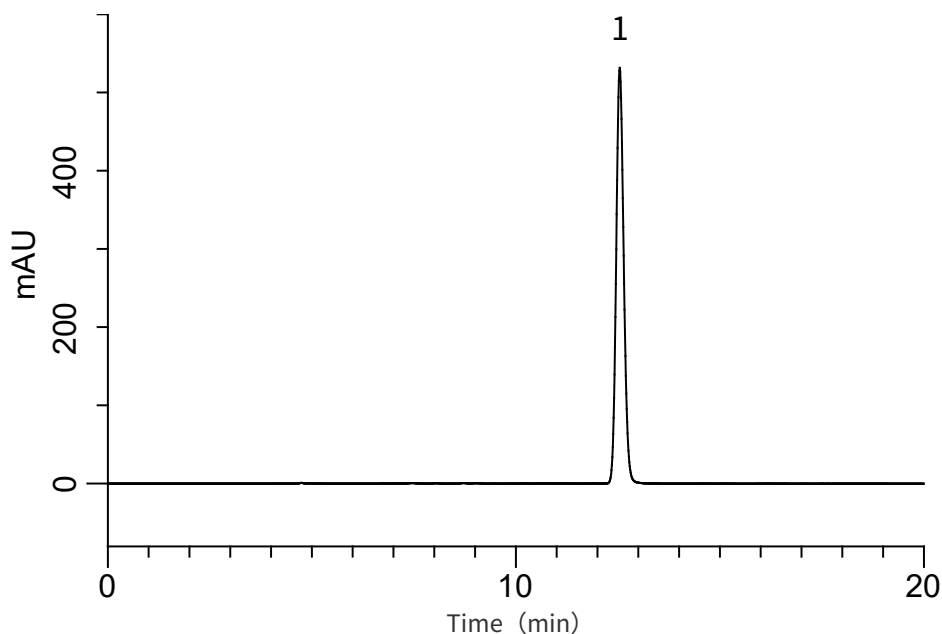

Conditions

System : GL7700 HPLC system
Column : InertSustainSwift C18
(5 μ m, 250 x 4.6 mm I.D.)
Column Cat. No. : 5020-88027
Eluent : A) CH₃OH
B) CH₃CN
C) 10 mM KH₂PO₄ in H₂O (pH 3.0, H₃PO₄)
A/B/C = 6/36/58, v/v/v
Flow Rate : 0.9 mL/min
Col. Temp. : 35 °C
Detection : UV 210 nm (PD7752 PDA Detector)
Injection Vol. : 20 μ L
Sample : Standard

Analyte:

1. Eplerenone 500 mg/L
Theoretical plates : 22,422 ($\geq$ 15,000)
Tailing factor : 1.12 ($\leq$ 1.5)
RSD of the
peak area (%) (n=6) : 0.06 ($\leq$ 1.0)