


Analysis of Tramadol Hydrochloride

(Under the Condition of the draft for the Japanese Pharmacopoeia)

Data No. LB472-0894

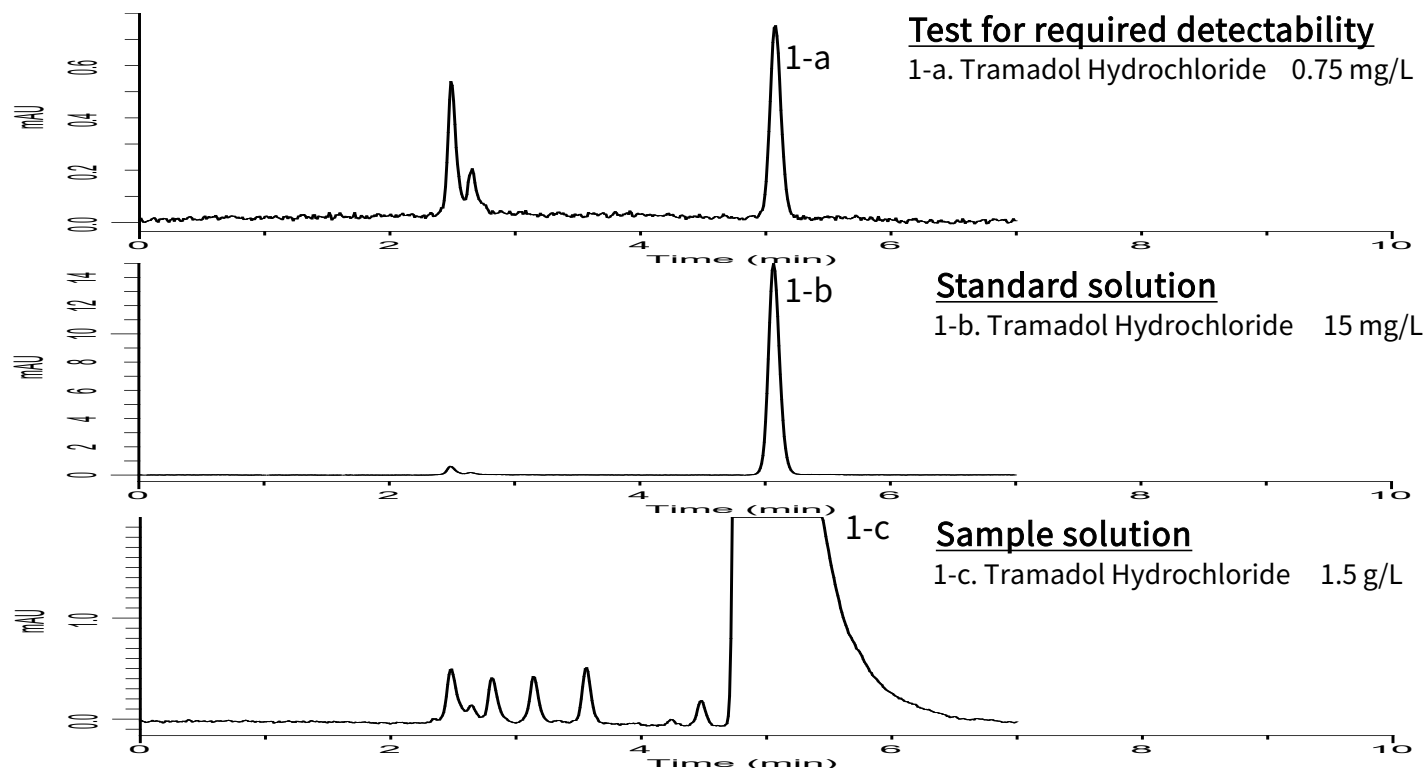

Conditions

System : GL7700 HPLC system
Column : InertSustainSwift C8
 (5 μ m, 250 x 4.0 mm I.D.)
Column Cat. No. : 5020-88324
Eluent : A) CH₃CN
 B) 0.2 % TFA in H₂O
 A/B = 59/141, v/v
Flow Rate : 0.9 mL/min
Col. Temp. : 25 °C
Detection : UV 270 nm (PD7752 PDA Detector)
Injection Vol. : 20 μ L
Sample : Standard

Analyte:

1. Tramadol Hydrochloride

The peak area ratio of 1-a to 1-b (%) : (3.5 $\leq$) 4.91 ($\leq$ 6.5)
 Theoretical plate number (1-b) : 15,016 ($\geq$ 5,000)
 RSD of the peak area (%) (n=6) : 0.20 ($\leq$ 2.0)
 Tailing factor (n=6) : 1.10 ($\leq$ 1.5)