


Analysis of Tulobuterol

(Under the Condition of the Japanese Pharmacopoeia 17th edition)

Data No. LB429-0894

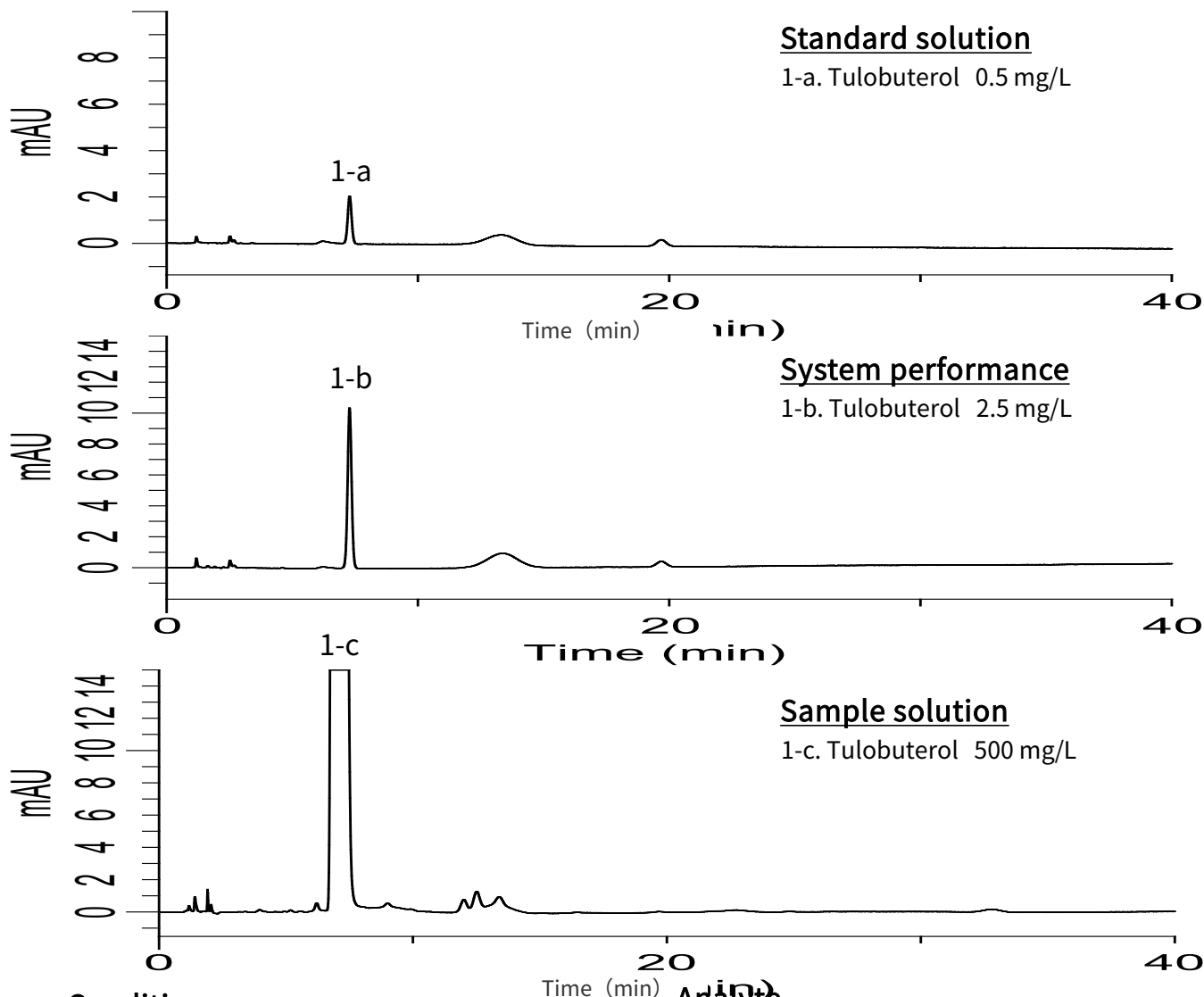

Conditions

System : GL7700 HPLC system
Column : InertSustain C18
(5 μ m, 150 x 4.6 mm I.D.)
Column Cat. No. : 5020-07345
Eluent : A) CH₃CN
B) Buffer*
A/B = 35/65, v/v

Flow Rate : 0.8 mL/min
Col. Temp. : 30 °C
Detection : UV 215 nm (PD7752 PDA Detector)
Injection Vol. : 25 μ L
Sample : Standard

*Dissolve 3.0 g of Sodium 1-octanesulfonate in 900 mL of water,
Add 5.0 mL of diluted phosphoric acid (1 $\rightarrow$ 150).

Analyte

1. Tulobuterol
Theoretical plate number : 10,026 ($\geq$ 5,000)
Tailing factor : 1.04 ($\leq$ 2.0)
RSD of the area (%) (n = 6) : 0.27 ($\leq$ 2.0)