


Analysis of Maltooligosaccharides

Data No. LB505-0874

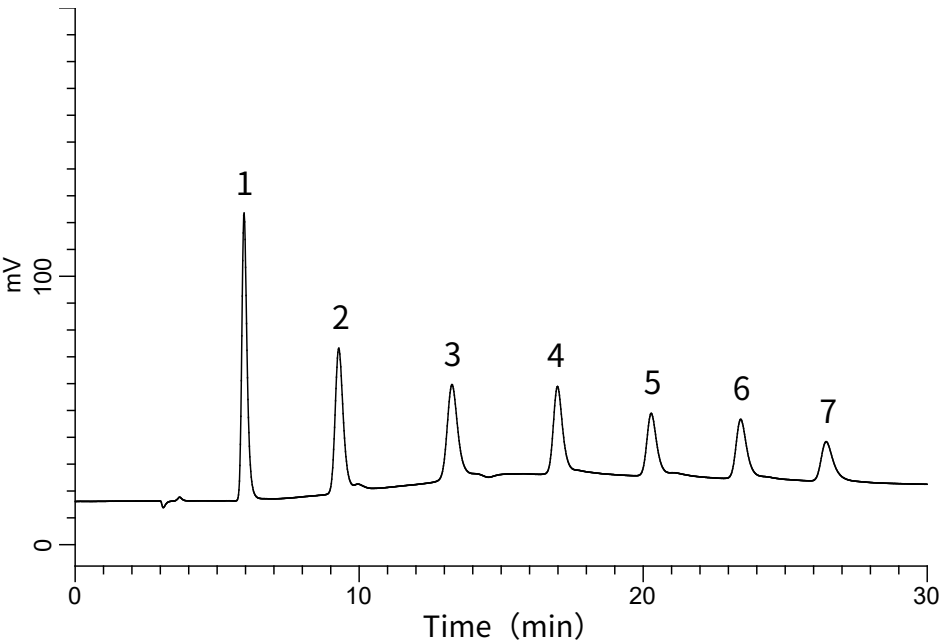

Conditions

System : GL7700 HPLC system
Column : InertSphere Sugar-1
(5 μm, 150 x 4.6 mm I.D.)
Column Cat. No. : 5020-11001
Eluent : A) 50 mM NaOH in H₂O
B) (50 mM NaOH + 400 mM CH₃COONa) in H₂O

Time (min)	A (vol%)	B (vol%)
0.0	85	15
30.0	50	50

Flow Rate : 0.5 mL/min
Col. Temp. : 30 °C
Detection : ECD Pulse Mode (ED723, Gold)
E1: 150 mV t1: 600 ms
E2: -1500 mV t2: 50 ms
E3: 600 mV t3: 50 ms
E4: -200 mV t4: 100 ms
ts: 50 ms

Injection Vol. : 10 μL
Sample : Standard

Analyte:

1. Glucose 10 mg/L
2. Maltose 10 mg/L
3. Maltotriose 10 mg/L
4. Maltotetraose 10 mg/L
5. Maltopentaose 10 mg/L
6. Maltohexaose 10 mg/L
7. Maltoheptaose 10 mg/L