


Cyromazine is triazine based molecule that is used as an insect growth inhibitor for the protection of livestock and poultry houses. Standard values have been set for its use in veterinary drugs, and a notification method using HPLC has been issued.


In this study, extraction and purification were performed with a mixed-mode polymer SPE (InertSep MCX FF) using a method based on the notification test method from the Ministry of Health, Labor and Welfare and literature examples\*, and analysis was performed using HPLC (detection at UV 240nm) and LC-MS/MS.

\*Tokyo Metropolitan Institute of Public Health Annual Report No.57 (2006)

## 1. SPE Pretreatment Flow

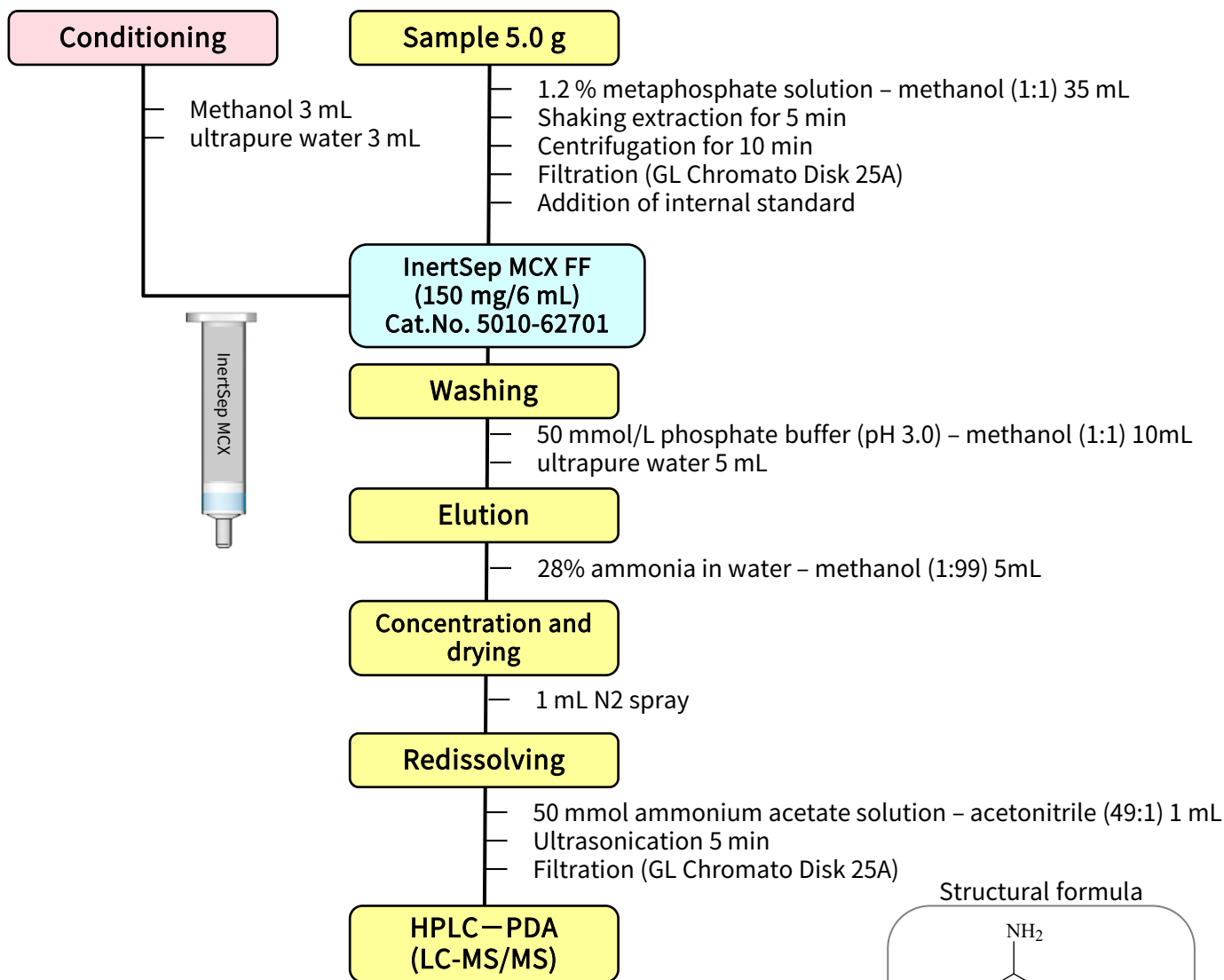


### 【InertSep MCX FF】

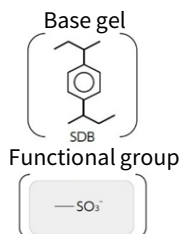

InertSep MCX FF is a styrene-divinylbenzene polymer (SDB) solid phase modified with a strong cation exchange group. By exerting both reverse phase and cation exchange actions, basic compounds can be strongly retained. Since the particle size is up to 70 μm, it is ideal for biological sample extracts.

Average Particle Size : 70 μm Pore Volume : 1.1 mL/g pH Range : 1-14  
Surface Area : 480 m<sup>2</sup>/g Pore diameter : 9nm

### Structural formula


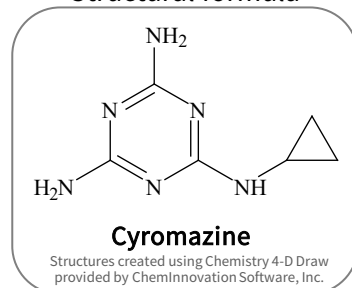


Recovery test with spiked extract (milk, 0.05 mg / kg, n = 2)

Rec (%)	HPLC-PDA	LC-MS/MS
Milk	80.1	78.6

## 2. Measurement conditions

### Standard Solution

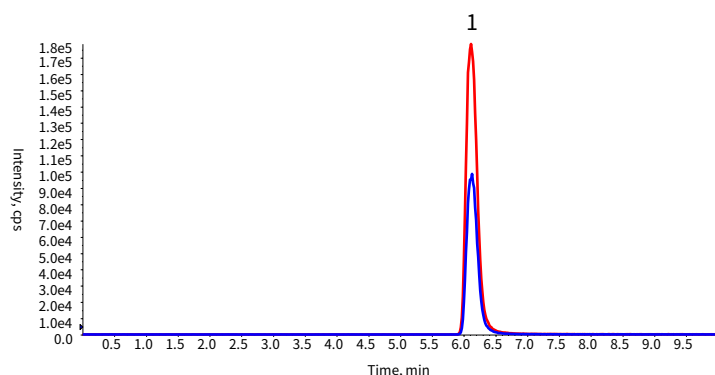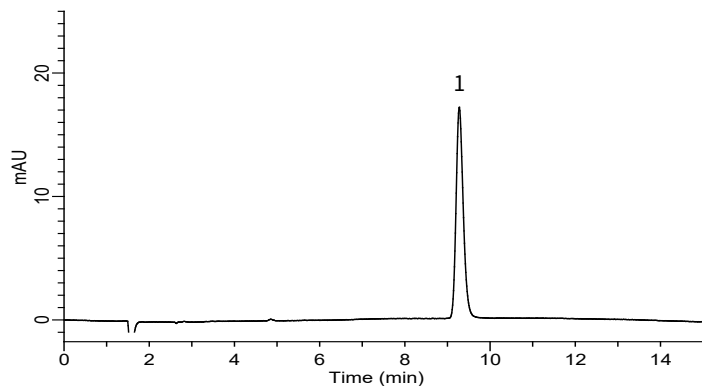

### HPLC conditions

#### <HPLC-PDA>

**Column** : InertSustain C18  
 (3 μm, 150 x 4.6 mm I.D.)  
**Column Cat. No.** : 5020-07445  
**Eluent** : A) 50 mmol/L CH<sub>3</sub>COONH<sub>4</sub>  
 B) CH<sub>3</sub>CN  
 A/B=49/1, v/v  
**Flow rate** : 1.0 mL/min  
**Col. Temp.** : 40 °C  
**Detection** : PDA 240 nm  
**Injection Vol.** : 50 μL  
**Sample** : 1. Cyromazine 1.75 mg/L

#### <LC-MS/MS>

**Column** : InertSustain C18  
 (3 μm, 150 x 2.1 mm I.D.)  
**Column Cat. No.** : 5020-14355  
**Eluent** : A) 50 mmol/L CH<sub>3</sub>COONH<sub>4</sub>  
 B) CH<sub>3</sub>CN  
 A/B=49/1, v/v  
**Flow rate** : 0.4 mL/min  
**Col. Temp.** : 40 °C  
**Detection** : LC/MS/MS  
 (4000QTRAP: ESI, Positive, SRM)  
**Injection Vol.** : 10 μL  
**Sample** : 1. Cyromazine 1.75 mg/L  
 Q1 / Q3 = 167.0 / 68.0 (Quantifier ion)  
 167.0 / 85.0 (Qualifier ion)

## 3. Applicability to other samples




Depending on the sample matrix, sufficient recovery of the target components may not be possible. If the matrix affects recovery, increasing the loading amount on the MCX cartridge may improve the recovery rate. However, if the loading amount is increased, peaks from interfering components will also increase, so it is important to select an appropriate loading amount suitable for the matrix.

### <Other samples: chicken>

InertSep MCX FF 150 mg/6 mL

InertSep MCX FF 250 mg/6 mL

InertSep MCX FF 500 mg/6 mL

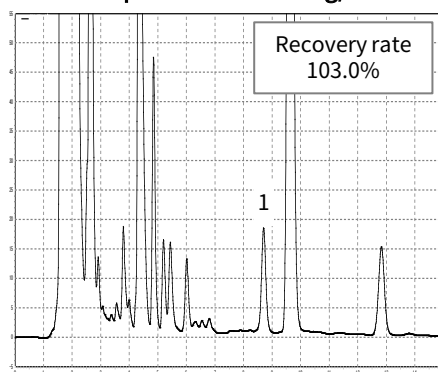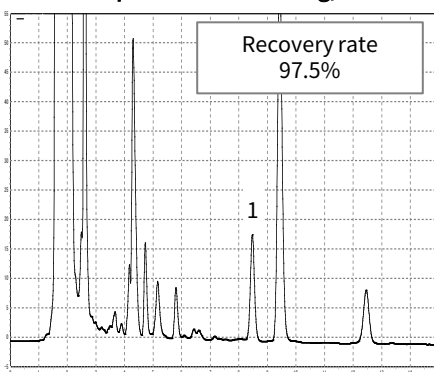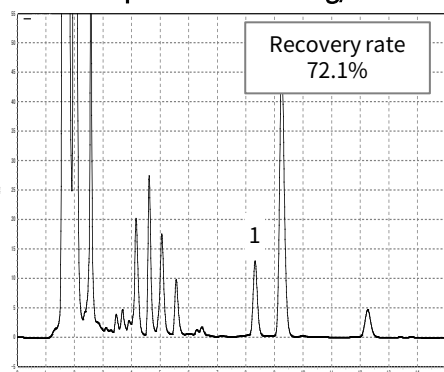

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

#### GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome  
Shinjuku-ku, Tokyo  
163-1130, Japan

Phone: +81-3-5323-6620  
Fax: +81-3-5323-6621  
Email: [world@glsc.co.jp](mailto:world@glsc.co.jp)  
Web: [www.glsciences.com](http://www.glsciences.com)

#### GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255  
Torrance, CA 90503  
USA

Phone: +1-310-265-4424  
Fax: +1-310-265-4425  
Email: [info@glsciencesinc.com](mailto:info@glsciencesinc.com)  
Web: [www.glsciencesinc.com](http://www.glsciencesinc.com)

#### GL Sciences B.V.

Dillenburgstraat 7C  
5652AM, Eindhoven  
The Netherlands

Phone: +31-40-254-9531  
Email: [info@glsciences.eu](mailto:info@glsciences.eu)  
Web: [www.glsciences.eu](http://www.glsciences.eu)

#### GL Sciences (Shanghai) Limited

Tower B, Room 2003  
Far East International Plaza  
No.317 Xianxia Road, Changning District  
Shanghai, China 200051

Phone: +86-21-62782272  
Email: [contact@glsciences.com.cn](mailto:contact@glsciences.com.cn)  
Web: [www.glsciences.com.cn](http://www.glsciences.com.cn)



International Distributors

Visit our Website at [www.glsciences.com/distributors](http://www.glsciences.com/distributors)